

**Epitomics RabMAb Supplement A
Product Data Sheet**

Catalog # EP401

Lot #: *Please refer to the vial*

Quantity: 50 ml

Biosafety Level: 1

Mycoplasma test*: Undetectable

Description –

RabMAb supplement A for hybridoma growth medium.

Protocols –

For complete growth medium:

Add 50 ml of Epitomics RabMAb Supplement A (containing antibiotic/antimycotic/Gentamicin) to 450 ml of RPMI 1640 medium with 0.05 mM 2-mercaptoethanol, and 10% fetal bovine serum. Incubate at 37°C.

Propagation:

Cultures can be propagated by centrifugation with subsequent re-suspension in fresh culture medium.

Establish new cultures at 1×10^5 cells/ml and maintain between 1×10^5 and 1×10^6 cells/ml.

Medium renewal: Every 2 to 3 days

Preservation: 90% FBS, 10% DMSO

Shipping/Storage Conditions: -20° C

Stability: 12 months from date of receipt

*Mycoprobe from R&D Systems (CUL001); VenorGeM from Sigma (MP0025)

For research use only. Not for use in diagnostic or therapeutic applications.

This product was manufactured under U.S. Patent No. 5,675,063. For a complete list of protocols and available related products, please visit www.epitomics.com.

Epitomics, Inc., 863 Mitten Road, Suite 103 Burlingame, California 94010-1303

Rev. A/0207/SL